Niskin Bottle NK

Manual de instrucciones / Instruction Manual

Botella Niskin NK

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
Niskin Bottle NK

The Niskin bottle is designed for the collection of water samples to the desired depth, activating its closure manually by messenger. It can be used individually or in series by attaching several bottles on the same line and actuating its closing by cascading messengers, allowing the collection of samples at various depths. The Niskin bottle is made of PVC, acetal, polyamide, stainless steel and latex. The tube contains no metallic elements inside. It is available in different volumes.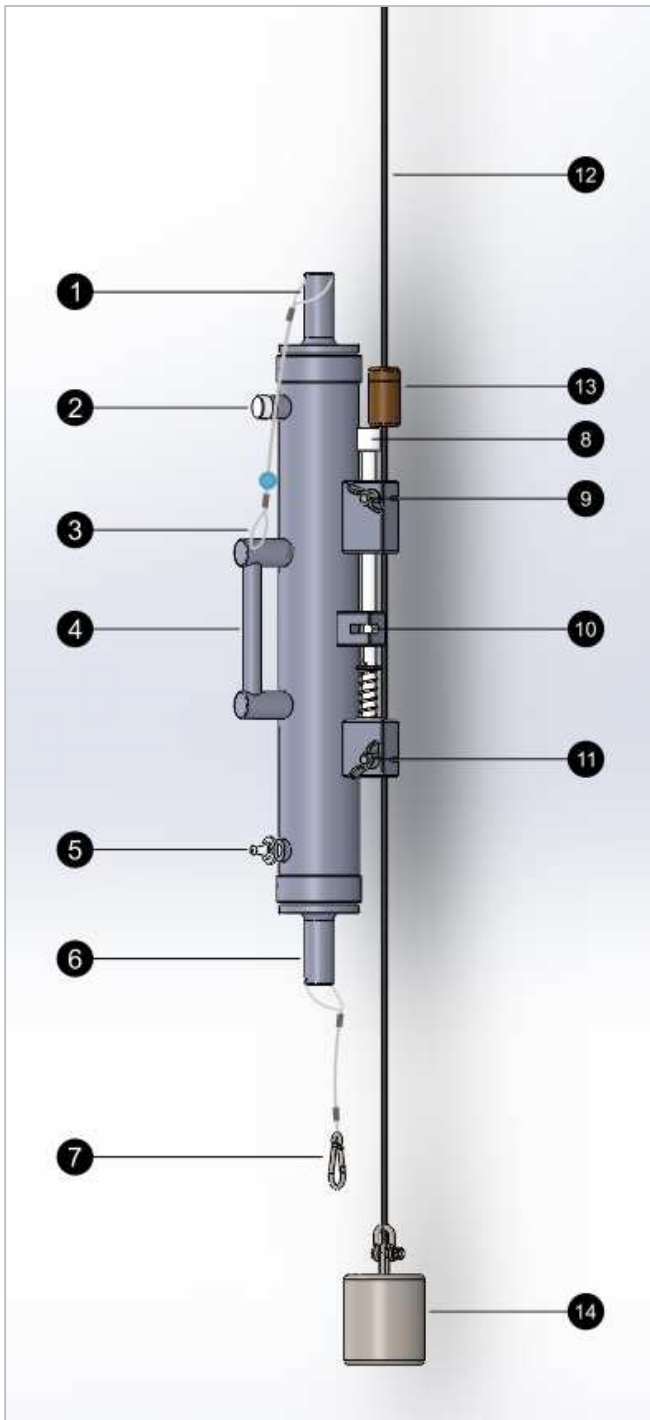


Niskin Bottle Components

1. Top plug.
2. Air screw.
3. Nylon with ball.
4. Handle.
5. Faucet.
6. Bottom plug.
7. Carabiner.
8. Impact base.
9. Upper fixing system.
10. Plugs releaser.
11. Lower fixing system.

Niskin Bottle Accessories

12. Main line.
13. Messenger.
14. Stainless steel cylindrical weight.

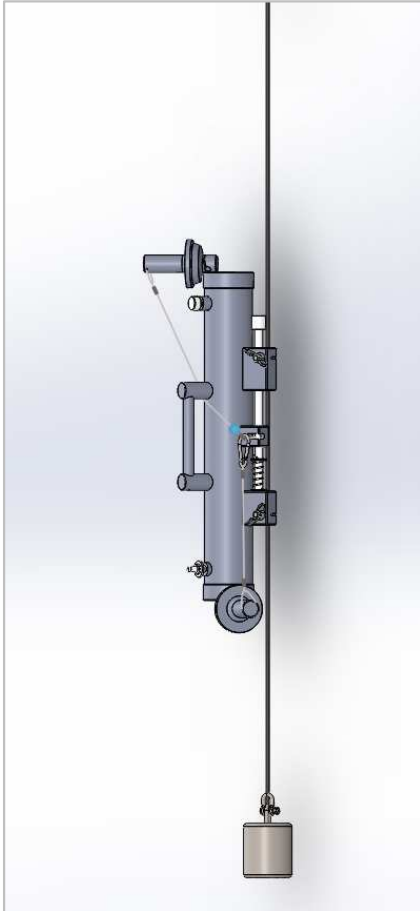Niskin Bottle fixing and loading

To fix the bottle to the main line (max. 8mmØ) loosen the fixing system (9 and 11) and place the line in the screw grooves. Tighten the wing nuts until firmly attaching the bottle to the line.

The bottle load is performed in the following steps and it is indicated on the bottle by adhesive vinyls:

- Pass the nylon with ball (3) through the handle (4), press the impact base (8) to open the plugs releaser (10) and insert the nylon into the metal pin (in case of using Niskin bottles in series also fix cascade messenger).
- Attach carabiner (7) to the nylon (3) between the ball and the plugs releaser (10).
- Ensure that the air screw (2) and the faucet (5) are closed. To close the faucet (5) pull it out.

Submerge the bottle to the desired depth.

Sample shooting and extraction

To close the bottle, place the messenger on the main line and drop it through it. The messenger will strike the impact base (8) and close the bottle.

Once closed, raise the bottle. To remove the sample, loosen the air screw (2) and open the faucet (5). The opening of the faucet (5) is performed by aligning the hole with the metal pin and pushing towards the bottle.

For more information contact:
info@aquaticbiotechnology.com